

#### DKG-21 M PERSONAL GAMMA RADIATION DOSIMETER

**Operating manual** BICT.412118.023-02 PЭ

#### CONTENTS

1 GENERAL GUIDELINES	5
2 MAIN INFORMATION ABOUT THE	
DOSIMETER	6
3 DESCRIPTION AND OPERATION	8
4 PROPER USE OF THE DOSIMETER	
5 TECHNICAL MAINTENANCE	65
6 STORAGE	

7 SHIPPING	102
/ SHIPPING	
8 DISPOSAL	
9 WARRANTY	
10 PACKING CERTIFICATE	
11 CERTIFICATE OF ACCEPTANCE	
12 REPAIR	111
13 ACCEPTANCE AND WARRANTY	
APPENDIX A	
APPENDIX B	115
APPENDIX C	116

This operating manual of BICT.412118.023-02 PЭ type is intended to inform the user about the principles of operation and rules of application of DKG-21 M personal gamma radiation dosimeter. The manual contains all information necessary for proper use of the DKG-21 M dosimeter (hereinafter called the dosimeter) and full realization of its technical possibilities.

Before operating the dosimeter, the user should be instructed on safety engineering and radiation safety, and should study this operating manual.

## **1 GENERAL GUIDELINES**

1.1 Carefully study the operating manual (hereinafter referred to as the OM) before using the dosimeter.

1.2 The OM should always be kept with the dosimeter.

1.3 All records in the OM should be accurate and clear. Notes made in pencil, erasures and uncertified corrections are not allowed.

# 2 MAIN INFORMATION ABOUT THE DOSIMETER

The dosimeter meets the TV V 33.2-22362867-010:2007 technical requirements.

The dosimeter is intended to be used as a standalone device or as part of the automated system of personal dosimetry control ASPDC-21 (hereinafter – the ASPDC) registered in the State Register for Measuring Instruments, and is accepted for application in Ukraine under No. Y1816-07.

Manufacturer:

Private Enterprise "SPPE "Sparing-Vist Center" 33 Volodymyr Velyky Str., Lviv 79026, Ukraine. tel.: (+38032) 242-15-15; fax: (+38032) 242-20-15. E-mail: sales@ ecotest.ua.

#### **3 DESCRIPTION AND OPERATION**

#### 3.1 Purpose of use

3.1.1 The dosimeter is designed to measure individual dose equivalent  $H_n(10)$  of gamma and Xray radiation (hereinafter - DE) and individual dose equivalent rate  $\dot{H}_{p}(10)$  (hereinafter – DER) of gamma and X-ray radiation, as well as support the automated database of the radiation burden to the personnel in the ASPDC

3.1.2 The dosimeter can be used at industrial enterprises and companies that deal with gamma sources.

#### **3.2 Technical specifications**

3.2.1 DER measurement range varies from 1.0  $\mu$ Sv/h to 1.0 Sv/h.

3.2.2 DER indication range of varies from  $0.01 \ \mu$ Sv/h to  $1.0 \ \mu$ Sv/h.

3.2.3 Main relative permissible error limit of DER measurement with 0.95 confidence probability is:

-  $\pm 20$  % - in the DER range of 1.0 to 10  $\mu$ Sv/h;

-  $\pm 15$  % - in the DER range of 10  $\mu$ Sv/h to 1.0 Sv/h.

3.2.4 DE measurement range varies from 0.001 to 9999 mSv.

3.2.5 Main relative permissible error limit of DE measurement with 0.95 confidence probability is  $\pm 15$  %.

3.2.6 Complementary relative permissible error limit of DER and DE measurements caused by ambient temperature change from -20 to +50 °C is  $\pm 5$  % per each 10 °C of deviation from +20 °C.

3.2.7 Energy range of registered gamma and X-ray radiation is from 0.05 to 6.00 MeV.

3.2.8 Energy dependence of DER and DE measurements in the energy range of 0.05 to 1.25 MeV is not more than  $\pm$  25 %.

3.2.9 Anisotropy at gamma quantum incidence at solid angle of  $\pm 60^{\circ}$  relative to the main measurement direction (perpendicular to the back panel of the dosimeter that is marked with the "+" symbol), not more than:

-  $\pm 25$  % - for <sup>137</sup>Cs and <sup>60</sup>Co radionuclides; -  $\pm 70$  % - for <sup>241</sup>Am radionuclide.

3.2.10 Time of operating mode setting during DER measurement is not more than 3 min.

3.2.11 Time of continuous operation of the dosimeter, which is charged from a new battery under conditions of normal radiation background, is not more than 1400 hours.

3.2.12 Unstable readings during an 8-hour continuous operation is not more than  $\pm 5$  %.

3.2.13 The dosimeter is powered from DC source of 2.4 to 3.2 V voltage and capacity of 560 mAh. Note – Nominal supply voltage – 3 V. 3.2.14 Current consumption of the dosimeter at nominal voltage under conditions of normal radiation background is not more than 0.4 mA.

3.2.15 Complementary relative permissible error limit of measurements caused by supply voltage deviation from the nominal voltage value, ranging from 3.2 to 2.4 V, is not more than  $\pm 10$  %.

3.2.16 Dimensions of the dosimeter is not more than  $98 \times 58 \times 18$  mm (without dimensions of a clip).

3.2.17 Weight is not more than 0.12 kg.

3.2.18 The dosimeter performs measurements under the following conditions:

- temperature – from - 20 to + 50  $^{\circ}$ C;

- relative humidity – up to 100 % at +  $25 \,^{\circ}$ C temperature;

- atmospheric pressure – from 60 to 106.7 kPa.

3.2.19 The dosimeter is resistant to the influence of atmospheric precipitation (rain).

3.2.20 The dosimeter is resistant to the influence of condensed atmospheric precipitation.

3.2.21 The dosimeter has "Clock" and "Alarm clock" operating modes.

3.2.22 The alarm clock rings during one minute.

3.2.23 DER and DE threshold level values with discreteness of a unit of a programmable digit over the whole operating measurement range can be programmed in the dosimeter from the personal computer (hereinafter – the PC) and in manual mode with the help of control buttons.

3.2.24 The dosimeter sends audio and light signals if the preset DER or DE threshold levels are exceeded.

3.2.25 DER and DE values as well as DER or DE threshold level values alternately appear on the digital liquid crystal display (hereinafter – the LCD), indicating the correspondence of information.

3.2.26 The history record of dose accumulation during the preset time interval is contained in the nonvolatile memory of the dosimeter with the discreteness of the record change of 1 min in the range of 5 to 255 min with reference to real time.

3.2.27 The dosimeter supports transmitting the history record of dose accumulation to the PC via the infrared port.

The distance of positive exchange between the dosimeter and the infrared port adapter (hereinafter – the IRPA) is in the range from 0.1 to 0.3 m.

3.2.28 With the help of the PC the dosimeter supports:

- blocking the mode of power supply switch off until data reading procedure is finished;

- blocking the modes of indication (DER, DER threshold, DE threshold, clock and alarm clock); change (DER threshold, DE threshold) and correction of clock and alarm clock settings.

3.2.29 The LCD can be switched off automatically not later than in 5 min after the switching-on of the dosimeter provided that the current gamma background is lower than the preset threshold level. At this point, none of the control buttons of the dosimeter should be pressed.

3.2.30 The dosimeter includes the mode of automatic power supply testing and its indication:

- power supply range from 2.5 to 2.6 V – segments (from 1 to 3) start blinking in the right upper corner of the LCD;

- power supply less than 2.4 V – four segments together with the indication of audio alarm on the LCD start blinking as well as audio alarm signal is annunciated.

3.2.31 Design of the dosimeter provides its decontamination.

3.2.32 Reliability requirements

3.2.32.1 Mean time to failure, not less than 10000 hours. Failure criterion is nonconformity to the requirements stated in 3.2.3.

3.2.32.2 Average operating life till the first major repair, not less than 10000 hours, and average service life, not less than 10 years.

Limit tolerance state criterion is deviation of parameters from the values stated in 3.2.3 that can not be eliminated.

3.2.32.3 Average shelf life is 10 years.

3.2.32.4 Average period of failure recovery at the authorized repair center or on the manufacturer premises without time required for shipping and verification after repair is not more than 2 hours.

#### **3.3 Delivery kit of the dosimeter** 3.3.1 The delivery kit is presented in Table 3.1. Table 3.1 – Delivery kit of the DKG-21 M dosimeter

Туре	Item	Quantity
BICT.412118.023-	Personal gamma	
02	radiation	1
	dosimeter	1
	DKG-21 M	
CR 2450	Battery	1

#### Table 3.1 (continued)

Туре	Item	Quantity
BICT.412118.023-	Operating	1
02 РЭ	manual	1
BICT.412915.018	Packaging	1
BICT.412914.004	Screwdriver	1
BICT.468353.014	USB/IrDA-	1 per delivery
	adapter	i per derivery
	CD Software	1 per delivery

# 3.4 Design and theory of operation

3.4.1 General information

The dosimeter is designed as a mono block with a built-in detector of gamma and X-ray radiation, with a printed circuit board equipped with a circuit of anode voltage formation, digital processing, control and indication, an infrared port of data exchange, and a battery.

The detector of gamma and X-ray radiation transforms gamma and X-ray radiation into the sequence of voltage pulses; the number of pulses is proportional to the registered radiation intensity. The circuit of anode voltage formation, digital processing, control and indication supports:

- scaling and linearization of the counting response of the detector;

- measurement of DER by measuring the average pulse frequency of the detector output;

- measurement of DE by measuring the total amount of pulses of the detector output;

- measurement of real time;

- formation and stabilization of anode voltage of the detector;

- control of operating modes of the dosimeter;
- indication of measurement results.

Power for operation is supplied by a disk-type lithium battery of CR 2450 (Panasonic).

#### 3.4.2 Design description

The dosimeter is designed as a flat rectangle plastic housing (Figure B), which consists of the front (1) and back (2) panels.

Three transparent windows are positioned on the upper part of the front panel. The LCD (3), the light-emitting diode indicator (4), as well as the optical system of the infrared port (5) are placed behind the windows.

Three acoustic holes (6) of the loudspeaker alongside the THRESHOLD (7) and MODE (8) control buttons are located in the middle of the front panel. The battery compartment, which is closed with the hermetic lid (9), is located at the bottom of the front panel. The clip (10) is on the back panel to fasten the dosimeter to clothes. The mechanical center (11) of the detector, being under the lid, is also inscribed there.

# 3.5 Labeling and sealing

3.5.1 Labeling is done in compliance with the requirements stated in design documents.

3.5.2 The front panel of the dosimeter is inscribed with the:

- full name of the device;

- trademark of the manufacturer;
- ingress protection rating of the housing IP54;
  approval mark.
- 3.5.3 The back panel of the dosimeter is inscribed with: - "Made in Ukraine";
- serial number of the dosimeter according to the numeration system of the producer enterprise;
  - mechanical center of the detector with "+" symbol;
  - year of manufacture.

3.5.4 Labeling quality is maintained during service life under all conditions and modes, except for labeling done on the individual package.

3.5.5 The dosimeter accepted by the Quality Control Department (QCD) and prepared for packing is sealed with a special film seal, covering the screw heads, which fasten together the upper and the lower panels, or a paste seal in the hollow above the head of the fastening screw.

## 3.6 Packing

3.6.1 Packing is performed in accordance with the requirements specified in design documents.

3.6.2 The dosimeter is packed in a special cardboard box, which (together with the operating manual), in its turn, is placed in a transparent polyethylene package, welded after packing performed.

# **4 PROPER USE OF THE DOSIMETER**

# 4.1 Operating limitations

4.1.1 Operating limitations are presented in Table 4.1.

Table 4.1 – Operating limitations

Name of operating limitations	Parameters of operating
limitations	limitations
1 Ambient air temperature	From - 20 to $+$ 50 °C
2 Relative humidity	Up to 100 % at 25 °C
2	temperature
3 Influence of gamma and	DER up to 10 Sv/h during 50 minutes
X-ray radiation	during 50 minutes

**4.2 Preparation of the dosimeter for operation** 4.2.1 Scope and order of external examination 4.2.1.1 Before using the dosimeter, unpack it and check if the delivery kit is complete. Examine for mechanical damages.

4.2.2 Rules and order of examination for operational readiness

4.2.2.1 Examine the control buttons before switching the dosimeter on.

4.2.2.2 Open the battery compartment and make sure the battery is inserted, connections are reliable, and there is no leakage of salts after durable storage of the dosimeter. In case there is salt leakage, remove the battery. Clean it, if possible, or replace, if not. Insert the battery and close the battery compartment.

4.2.3 Guidelines on switching on and testing the dosimeter

4.2.3.1 Prepare the dosimeter for operation. Do the following:

- unpack the dosimeter;

- open the battery compartment and insert the battery of CR 2450 type observing the polarity. The dosimeter should enter the mode of DER measurement indicated by " $\mu$ Sv/h" measurement units, which are constantly displayed on the LCD.

4.2.3.2 Press shortly the MODE button and make sure the dosimeter has entered the mode of DE indication. DE units of measurement expressed in "mSv" should appear on the LCD. 4.2.3.3 Press shortly the MODE button and make sure the dosimeter has entered the mode of real time indication, which is shown by a one-second blinking colon between the two pairs of the LCD digits.

4.2.3.4 Press shortly the MODE button and make sure the dosimeter has entered the mode of the alarm clock setting, which is indicated by an unblinking colon between the two pairs of digits on the LCD. After setting the alarm clock (4.3.3.7) press shortly the MODE button to switch it on, which should be indicated by a ")))" symbol on the LCD. 4.2.3.5 Hold the MODE button pressed for 4 s to switch the dosimeter off.

**Note.** The battery should be replaced if the low battery symbol appears on the LCD – gradual blinking of one to four segments of the battery symbol in the right top corner of the LCD, which depends on the increasing process of discharge, when the dosimeter is switched on irrespective of the selected mode. All four segments of the symbol are blinking and a short audio signal is generated each 4 s when the battery is fully discharged.

4.2.4 **List of possible troubles and troubleshooting** 4.2.4.1 The list of possible troubles and troubleshooting is presented in Table 4.2.

Table 4.2 - Tossible troubles and troubleshooting		
Trouble	Probable cause	Troubleshooting
1 The	1 The battery is	1 Replace the
dosimeter	discharged	battery
does not	2 Poor contact	2 Rest ore the
switch on at	between the battery	contact between
pressing the	and the battery	the battery and the
MODE button	compartment clamps	clamps

Table 4.2 - Possible troubles and troubleshooting

## Table 4.2 (continued)

Trouble	Probable cause	Troubleshooting
2 The "Err" symbol is displayed on the LCD after the battery has been replaced	Failure of the nonvolatile memory of the dosimeter	Send the dosimeter for repair to the manufacturer

Table 4.2 (continued)

Trouble	Probable cause	Troubleshooting
3 The "Err1"	Failure of the	Send the
symbol is	anode voltage	dosimeter for
displayed on the	former or the	repair to the
LCD during	ionizing	manufacturer
operation of the	radiation	
dosimeter	detector	

4.2.4.2 At failure to eliminate the troubles presented in Table 4.1, or at detection of more complicated troubles, the dosimeter should be sent for repair to the manufacturer.

### 4.3 Use of the dosimeter

4.3.1 Safety measures

4.3.1.1 The dosimeter meets the requirements of the valid standards with regard to people protection against electric-shock hazard.

4.3.1.2 The dosimeter contains electric circuits of voltage up to 400 V; demount the dosimeter when the power supply is switched off.

4.3.1.3 The dosimeter's design excludes any electric voltages on the outside of the dosimeter.

4.3.1.4 A special protective jacket is used to prevent accidental contact with conductive parts.

4.3.1.5 Ingress protection rating is IP54.

4.3.1.6 Disposal of the dosimeter should be performed in compliance with the general rules, i.e. metal is recycled or melted, and plastic parts are dumped.

**Note.** If the dosimeter is contaminated by any liquid or dry radionuclides and it is impossible to completely deactivate the device, the dosimeter should be buried as solid radioactive waste. 4.3.2 Operating modes of the dosimeter

4.3.2.1 The dosimeter operates in the following modes:

- switching the dosimeter on/off;
- measurement and indication of DER;
- programming of audio and light alarm threshold levels of DER;
  - indication of DE to be measured;

- programming of audio and light alarm threshold levels of DE;

- indication and correction of real time;

- indication and correction of the alarm clock settings, switching the alarm clock on/off;

- power supply control;

- performance monitoring of the ionizing radiation detector.

#### 4.3.3 Operation procedure of the dosimeter 4.3.3.1 Switching the dosimeter on/off

Press shortly the MODE button to switch the dosimeter on. The information displayed on the LCD shows that the dosimeter is on.

At the same time the infrared port of the dosimeter is activated for 15 s. While the infrared port is active, you can perform the procedure of data exchange with the PC, programming of the threshold levels, intervals of dose history accumulation, and permit/prohibit certain operating modes of the dosimeter. The blinking digits on the LCD indicate that the infrared port is active. As soon as data exchange with the PC is finished, the dosimeter starts accumulating dose history with the preset interval. Otherwise, the dosimeter starts operating

in the stand-alone mode with integral dose accumulation without dose history accumulation.

Press the MODE button once again and hold it pressed for 4 s to switch the dosimeter off. If the dosimeter is switched on in the stand-alone mode, e.i. no data exchange with the PC is done, the dosimeter will be switched off. If during switching it on data exchange between the dosimeter and the PC took place the effort to switch the dosimeter off would only activate the infrared port for 15 s.

### 4.3.3.2 Measurement of DER

The mode of DER measurement is entered automatically after the dosimeter is switched on. The units of measurement are expressed in µSv/h. The process of DER measurement accumulation and averaging will start after the dosimeter is switched on. The process will continue up to 1600 s at DER values close to background. The data on the LCD will be updated each 10 s. However, close to true result will appear in 2-3 minutes. At increase of the radiation intensity, the time of DER measurement averaging and the time of data updating on the LCD will decrease to minimum 2 s.

The units of measurement are expressed in  $\mu$ Sv/h, mSv/h, and Sv/h.

The statistical error of the displayed DER measurement result is indicated by the blinking or steady decimal point. The blinking point indicates that the statistical error of the indicated DER measurement result exceeds maximum permissible error. Hence, the measurement result can be used only for rough DER evaluation. The steady point informs that the statistical error of the DER measurement result is within the permissible range.

The main direction of the dosimeter when DER is measured is the direction perpendicular to the front (rear) panel of the dosimeter.

DER measurement result is considered to be the arithmetic mean of five last measurements in 8 min after the intensity of radiation is changed at DER levels in the range of 1.0 to 10.0  $\mu$ Sv/h, or within 2 min to 2 s for the levels in the range of 10.0  $\mu$ Sv/h to 1.0 Sv/h.

Measurement intervals and subranges will be set automatically according to the intensity of registered radiation. **Note.** The process of data averaging can be stopped forcefully for quick estimating of DER. To do this, press shortly the THRESHOLD button. Rough evaluation of DER value will be shown within 1 min.

## 4.3.3.3 **Programming of audio and light DER alarm threshold levels**

Audio and light alarm threshold levels of DER are programmed in the mode of DER measurement. Press the THRESHOLD button and hold it pressed for 4 s to start programming. The low-order digit will start blinking on the LCD. Set an appropriate value of the low-order digit by pressing and releasing the THRESHOLD button. Press shortly the MODE button to proceed to the programming of another digit, the latter will start blinking. The necessary value of the digit is set with short pressing and releasing the THRESHOLD button.

After the last digit of the threshold level is set and at next pressing of the MODE button, the LCD will blink four times, indicating that the value is fixed. The dosimeter will enter the mode of DER measurement. Press the THRESHOLD button and hold it pressed not more than 4 s after a new threshold level value appears to check the value of the fixed DER threshold level.

When holding the THRESHOLD button pressed for more than 4 s the low-order digit starts blinking, indicating that a new threshold level value can be programmed.

Blinking of the red light-emitting diode and a twotone audio alarm indicate that the programmed DER threshold level has been exceeded.

### Notes

1 The DER threshold level value of 1.0  $\mu$ Sv/h is set automatically after the dosimeter is switched on.

2 A preset zero value of the DER threshold level sets off the alarm system when the threshold level is exceeded.

### 4.3.3.4 Indication of DE measurement

Press shortly the MODE button to enter the mode of DE measurement indication. This mode follows the mode of DER measurement (entered automatically after the dosimeter is switched on).

The "mSv" symbol that appears on the LCD indicates you have entered the appropriate mode. A comma after the first left digit will appear on the LCD after the dosimeter is switched on The comma will automatically shift to the right until full completion of the DE scale of the dosimeter as the DE measurement value increases. The units of DE measurement are expressed in mSv.

# 4.3.3.5 Programming of audio and light DE alarm threshold levels

Audio and light alarm threshold level of DE is programmed in the mode of DE measurement. Hold the THRESHOLD button down for more than 4 s to start programming. The low-order digit will start blinking on the LCD.

Set an appropriate value of the low-order digit by pressing and releasing the THRESHOLD button. Press shortly the MODE button to proceed to programming of another digit, the latter will start blinking. Set an appropriate value of the digit by pressing and releasing the THRESHOLD button.

After the last digit of the threshold level is set and at next pressing of the MODE button, the LCD will blink four times, indicating that the value is fixed, and the dosimeter will enter the mode of DE indication.

Press the THRESHOLD button and hold it pressed for maximum 4 s after the threshold level value appears to check the value of the fixed DE threshold level. Hold the THRESHOLD button pressed for more than 4 s to set the threshold value to zero. The low-order digit will start blinking at that indicating that a new threshold level value can be programmed.

Blinking of the red light-emitting diode and a twotone audio alarm indicate that the preset DE threshold level has been exceeded.

At reaching 90% of the programmed threshold level, the dosimeter sends an interrupted audio signal to inform the user about possible near reaching the DE threshold level. Press any button to switch this audio alarm off. **Note.** The DE threshold level value of 0.000 mSv is set automatically when the dosimeter is switched on indicating that the alarm is switched off.

## 4.3.3.6 Indication and correction of real time

Press shortly the MODE button to enter the mode of real time indication. This mode follows the mode of indication of DE measurement.

It is indicated by a one-second blinking ":" symbol between the two pairs of the LCD digits.

The digits from the right to the left show the following: the first digit indicates minutes; the second one - tens of minutes; the third one - hours; the fourth one - tens of hours.

Press the THRESHOLD button and hold it down until two digits to the right from the ":" symbol start blinking to correct the value of real time. Release the button afterwards. The proper values of units and tens of minutes are fixed by further pressing and holding the THRESHOLD button The minutes can also be corrected by short pressing of the THRESHOLD button. Each pressing will change the value per unit. Press shortly the

MODE button to correct the value of hours. The two digits to the left from the ":" symbol start blinking at that. The hour value correction is performed likewise. Press shortly the MODE button once again to exit the mode of real time correction.

## 4.3.3.7 Indication and correction of alarm clock settings, switching alarm clock on/off

Press shortly the MODE button to enter the mode of indication of the alarm clock time. This mode follows the mode of real time indication. A non-blinking ":" symbol between the two pairs of digits on the LCD indicates you have entered the appropriate mode.

Press the THRESHOLD button and hold it down until the two digits to the right from the ":" symbol start blinking to correct the alarm clock settings. Release the button afterwards. Set the proper values of units and tens of minutes by further pressing and holding the THRESHOLD button The minutes can also be corrected by short pressing of the THRESHOLD button. The value will change per unit each time in this case. Press shortly the MODE button to correct the value of hours. The two digits to the left from the ":" symbol start blinking at that. Hour values can be corrected likewise.

Press shortly the MODE button to switch the alarm clock on/off after correcting the hour values of the alarm clock. A blinking sound symbol ")))" should appear on the LCD. Press shortly the THRESHOLD button; make the non blinking sound symbol appear on the LCD to switch the alarm clock on.

Press shortly the THRESHOLD button to switch the alarm clock off. The sound symbol should extinct. Fix the alarm clock settings by further short pressing of the MODE button. If the alarm clock is on, the sound symbol will be displayed on the LCD irrespective of the selected operating mode.

Note. The alarm clock will continue working even after the power supply of the dosimeter is off (provided that the batteries are inserted). The dosimeter will automatically enter the mode of real time indication when the alarm clock goes off. Press any control button to switch off audio signal of the alarm clock. Otherwise, audio signal will be disabled automatically in a minute after the alarm clock rings.

### 4.3.3.8 Power supply control

The mode of power supply control is turned on automatically at switching on the dosimeter. It is

indicated by the four-segment battery symbol displayed in the right upper corner of the LCD. The number of segments blinking starting from the last right one shows the level of battery discharge. Blinking of three and more segments means that the battery should be replaced with new one.

### 4.3.3.9 Operability check of the detector

The mode of detector's operability check is switched on as soon as the dosimeter is on. If the detector fails, the "Err1" symbol is displayed on the LCD, which means the dosimeter should be sent for repair.

### **5 TECHNICAL MAINTENANCE**

## 5.1 Technical maintenance of the dosimeter

5.1.1 General instructions

The list of operations performed during technical maintenance (hereinafter called TM) of the dosimeter, the order and the peculiarities of operational phases are presented in Table 5.1.

# Table 5.1 - List of operations during technical maintenance

		TM type		ſ
	during		During	OM item No.
Operations	everyday	periodical	long-	Νο
	use	use	term	NO
	use	(annually)	storage	Ũ
External examination	+	+	+	5.1.3.1
Delivery kit completeness check	-	+	+	5.1.3.2
Operability check	+	+	+	5.1.3.3
Power supply disconnection	-	+	+	5.1.3.4
Verification of the dosimeter	-	+	+	5.2
Note. "+" means the operation is applicable for this type of TM;				
"-" means the operation is not applicable.				

5.1.2 Safety measures

5.1.2.1 TM safety measures fully comply with safety measures stated in item 4.3.1 of the present OM.

5.1.3 Maintenance procedure of the dosimeter

5.1.3.1 External examination

External examination of the dosimeter should be performed in the following order:

- check the technical condition of surface, inspect for integrity of seals, absence of scratches, traces of corrosion, surface damage of the dosimeter;

- check the condition of clamps in the battery compartment.

5.1.3.2 Delivery kit completeness check

Check if the delivery kit is complete according to Table 3.1.

5.1.3.3 Operability check of the dosimeter.

Operability check of the dosimeter is performed according to item 4.2.3 of the present OM.

5.1.3.4 Power supply switch off

Power supply should be switched off before the long-term storage of the dosimeter. Do this as follows:

- switch the dosimeter off;
- open the lid of the battery compartment;
- remove the batteries;

- examine the battery compartment, check the contact clamps accuracy, clean the battery compartment from contamination and contact clamps from oxides;

- make sure there is no humidity, no salt spots on the surface of the batteries, and no damages of the insulated coating.

### 5.2 Verification

5.2.1 The DKG-21 M dosimeter should be verified after manufacture, repair or during use.

5.2.2 The interval between verifications should not exceed 12 months.

5.2.3 Verification operations are presented in Table 5.2.

## Table 5.2 - Verification operations

Operation name	Verification technique No.
External examination	5.2.7.1
Testing	5.2.7.2
Calculation of main relative permissible error limit of DER measurement in the DER range of $1.0 \mu$ Sv/h to $1.0 $ Sv/h	5.2.7.3, 5.2.7.4
Calculation of main relative permissible error limit of DE measurement	5.2.7.3, 5.2.7.5
Presentation of verification results	5.2.7.6

5.2.4 Verification facilities are presented in Table 5.3. Table 5.3 - Verification facilities

Name	Regulatory Document or Main Technical Specifications
УПГД-3Б	DER range from 0.01 µSv/h to 1 Sv/h. Energy range from 59 KeV to 1.25 MeV.
testing	
equipment	Main relative permissible error limit of DER
	and DE measurements is 4 % at 0.95
	confidence probability

### Table 5.3 (continued)

Name	Regulatory Document or Main Technical Specifications
Phantom	Dimensions: $30 \times 30 \times 15$ cm; PMMA walls (polymethylmethacrylate, front wall thickness – 2.5 mm, other walls thickness – 10 mm); phantom is filled with distilled water
Aspiration psychrometer	Temperature measurement range from - 30 °C to + 50 °C. Temperature measurement error is $\pm 0.1$ °C. Relative humidity measurement range from 10 to 100 %. Relative humidity measurement error from $\pm 12$ % at t = -10 °C to $\pm 2$ % at t = 30 °C

Table 5.3 (continued)

ruble 5.5 (continued)			
Name	Regulatory Document or Main Technical Specifications		
Stop-watch	Measurement range from 1 s to 59 min		
Control aneroid barometer	Pressure measurement range from 81.3 to 105.3 kPa (from 610 to 790 mm Hg). Pressure measurement error is $\pm 0.107$ kPa (0.8 mm Hg)		
Notes 1 All verification facilities should be certified, tested or calibrated. 2 Use of other measuring instruments, tools and equipment with specifications similar to those outlined in Table 5.3 is allowed.			

5.2.5 Verification should be performed in accordance with safety measures presented in 4.3.1 of the OM.

5.2.6 Verification conditions

Verification should be performed under the following conditions:

- ambient air temperature range within (20±5) °C;
- relative air humidity from 30 to 80 %;
- atmospheric pressure from 86 kPa to 106.7 kPa;

- natural background level of gamma radiation should not exceed  $0.3 \ \mu$ Sv/h;

- power supply voltage within  $(3.0\pm0.2)$  V.

# 5.2.7 Verification procedure

5.2.7.1 External examination

5.2.7.1.1 During external examination the dosimeter should meet the following requirements:

- the delivery kit should be completed as stated in Table 3.1;

- labeling should be accurate;
- QCD seals should not be violated;

- the dosimeter should be free from mechanical damage that may affect its performance.

5.2.7.1.1.2 If the requirements in 5.2.7.1.1 are satisfied, proceed to the next verification operation.

5.2.7.1.1.3 If the delivery kit is not completed as stated in Table 3.1, verification should be stopped until the delivery kit is complete.

5.2.7.1.1.4 If labeling and sealing requirements are not satisfied, and the dosimeter has the signs of mechanical damages that affect its performance, it can not be verified and should be sent for repair. 5.2.7.2 Testing

5.2.7.2.1 Perform operations stated in 4.2.3.

5.2.7.2.1.1 If all operations stated in 4.2.3 are performed, proceed to the next test operation.

5.2.7.2.1.2 Even if a single operation stated in 4.2.3 can not be performed, the dosimeter should not be verified and should be sent for repair.

5.2.7.3 DER and DE measurements should be performed on the phantom with 30x30x15cm dimensions, with PMMA walls (polymethylmethacrylate, front wall thickness - 2.5 mm, other walls thickness - 10 mm); the phantom should be filled with distilled water.

5.2.7.3.1 During measurement the dosimeter should be placed close to the phantom surface, directed to gamma source. The indicator of the dosimeter should be directed towards gamma source.

5.2.7.4 Calculation of main relative permissible error limit of DER measurement in the DER range of  $1.0 \,\mu$ Sv/h to  $1.0 \,$ Sv/h is performed as follows.

5.2.7.4.1 Prepare the dosimeter for DER measurement and program DER threshold level to the value of 1  $\mu$ Sv/h.

5.2.7.4.2 Fix the dosimeter on the phantom and the УПГД-3Б carriage so that the mechanical center of the УПГД-3Б collimator coincides with the mechanical center of the dosimeter's detector that is marked on the back panel of the dosimeter. Wait for 8 s and proceed to take ten measurements of background DER ( $\dot{H}_{ndi}$  (10))

in УПГД-3Б with 10 s interval. Calculate the average DER value in  $\mu$ Sv/h by the formula:

$$\overline{\dot{H}}_{p\phi}(10) = \frac{\sum_{i=1}^{10} \dot{H}_{p\phi i}}{10}$$
(5.1)

5.2.7.4.3 Place the УΠΓД-3Б carriage together with the phantom and the dosimeter in the position, where DER from the <sup>137</sup>Cs source with radionuclide is  $\dot{H}_{p0}(10) = (1.15 \pm 0.15) \,\mu$ Sv/h. After irradiation of the dosimeter was started, wait for 3 min and proceed to take ten measurements of DER with 10 s interval. Calculate the average DER value  $(\overline{H}_{p\Sigma}(10))$  by the formula (5.1). Calculate the DER value without gamma background DER of the УПГД-3Б in  $\mu$ Sv/h by the formula

$$\overline{\dot{H}}_{p}(10) = \overline{\dot{H}}_{p\Sigma}(10) - \overline{\dot{H}}_{p\phi}(10)$$
(5.2)

Note - The distance between the mechanical center of the source and the mechanical center of the dosimeter's detector is considered to be the distance between the mechanical center of the source and 81 the plane, which is perpendicular to the direction of gamma-quanta beam propagation, and passes through the mechanical center of the dosimeter in this plane.

5.2.7.4.4 Check if the dosimeter sends audio and light signal.

5.2.7.4.5 Calculate main relative permissible error limit of DER measurement in percentage by the following formula:

$$\Delta = 1, 1 \sqrt{\delta \dot{H}_{p}(10)^{2} + \delta \dot{H}_{p0}(10)^{2}}, \qquad (5.3)$$

where  $\delta \dot{H}_p(10)$  – relative error of measurement result calculated by the formula:

$$\delta \dot{H}_{p}(10) = \frac{\dot{H}_{p}(10) - \dot{H}_{p0}(10)}{\dot{H}_{p0}(10)}, \qquad (5.4)$$

where  $\delta \dot{H}_0(10)$  – main relative permissible error limit of DER measurement from the gamma reference source.

5.2.7.4.6 Set the DER threshold level value to 10  $\mu$ Sv/h and perform operations 5.2.7.4.3-5.2.7.4.5 for DER  $\dot{H}_{p0}(10) = (12 \pm 2) \mu$ Sv/h in 1 min after the dosimeter is exposed to radiation provided that DER measurement time is 5 s, n = 5.

5.2.7.4.7 Set the DER threshold level value to 100  $\mu$ Sv/h and perform operations 5.2.7.4.6 for DER  $\dot{H}_{p0}(10) = (120 \pm 20) \,\mu$ Sv/h.

5.2.7.4.8 Set the DER threshold level value to 1 mSv/h and perform operations 5.2.7.4.6 for DER  $\dot{H}_{p0}(10) = (1.2 \pm 0.2) \text{ mSv/h}.$ 

5.2.7.4.9 Set the DER threshold level value to 10 mSv/h and perform operations 5.2.7.4.6 for DER  $\dot{H}_{p0}(10) = (12 \pm 2) \text{ mSv/h}.$ 

5.2.7.4.10 Program DER threshold level to the value of 100 mSv/h and perform operations 5.2.7.4.6 for DER  $\dot{H}_{p0}(10) = (120 \pm 20)$  mSv/h.

5.2.7.4.11 Set the DER threshold level value to 800 mSv/h and perform operations 5.2.7.4.6 for DER  $\dot{H}_{p0}(10) = (900 \pm 100) \text{ mSv/h}.$ 

5.2.7.4.12 Maximum value of all received errors is the limit of main relative permissible error in DER measurement.

5.2.7.4.13 If the main relative permissible error limit of DER measurement with 0.95 confidence probability does not exceed:

-  $\pm 20$  % in the DER range of 1.0 to 10  $\mu$ Sv/h;

-  $\pm 15$  % in the DER range of 10  $\mu$ Sv/h to 1.0 Sv/h, proceed to the following verification operation.

5.2.7.4.14 If the main relative permissible error limit of DER measurement does not meet the requirements stated in 5.2.7.4.13, the dosimeter can not be verified and should be sent for repair.

5.2.7.5 Calculation of main relative permissible error limit of DE measurement is done in the following way.

5.2.7.5.1 Switch the dosimeter on and set the DER threshold level value to  $0.00 \,\mu$ Sv/h, the DE threshold level value –  $0.001 \,\text{mSv/h}$ . Prepare the dosimeter for DE measurement. The initial DE readings should be  $0.000 \,\text{mSv}$ .

5.2.7.5.2 Fix the dosimeter on the phantom and the  $\Psi\Pi\Gamma\Pi_{-35}$  carriage so that the mechanical center of the  $\Psi\Pi\Gamma\Pi_{-35}$  collimator coincides with the mechanical center of the dosimeter's detector that is marked on the back panel of the dosimeter.

5.2.7.5.3 Place the УΠΓД-3Б carriage together with the phantom and the dosimeter in the position, where DER from the <sup>137</sup>Cs source with radionuclide is  $\dot{H}_{p0}(10) = (1.15 \pm 0.15) \,\mu$ Sv/h and at the same time switch on the standard stop-watch. Proceed to direct the source towards the collimator.

5.2.7.5.4 In the period of 7200 s measurement time take results of DE. Check if audio and light alarms were actuated by this. After that switch the dosimeter off.

5.2.7.5.5 Switch the dosimeter on and set the threshold level value of DER to  $0.00 \,\mu\text{Sv/h}$ , the threshold level value of DE  $- 0.01 \,\text{mSv/h}$ .

5.2.7.5.6 Prepare the dosimeter for DE measurement. Place the  $Y\Pi\Gamma\Pi$ -35 carriage together with the phantom and the dosimeter in the position, where DER from the <sup>137</sup>Cs source with radionuclide is  $\dot{H}_{p0}(10) = (12 \pm 2) \mu$ Sv/h and at the same time switch on the standard stop-watch. Proceed to direct the source towards the collimator. 5.2.7.5.7 In the period of measurement time calculated by the formula  $t = 3600 + t_{\partial}$ , c, where  $t\partial$ , c is the period of time when the source is directed towards the collimator, take DE measurement results. Check if the dosimeter sends audio and light signals. Switch the dosimeter off afterwards.

5.2.7.5.8 Calculate main relative permissible error limit of DE measurement in percentage by the formula

$$\delta H_p(10) = 1,1 \sqrt{\left(\frac{H_p(10) - H_{p0}(10)}{H_{p0}(10)}\right)^2 + \left(\frac{\delta H_{p0}(10)}{2}\right)^2}, \quad (5.5)$$
  
where  $H_{p0}(10) = \dot{H}_{p0}(10) \cdot t$  – DE of the VIIITД-36 carriage;  
 $\delta H_{p0}(10)$  - main relative permissible error limit of DE measurement of the VIIITД-36 carriage is defined as follows:

$$\delta H_{p0}(10) = \sqrt{\left(\delta \dot{H}_{p0}(10)\right)^2 + \left(\delta t\right)^2} , \qquad (5.6)$$

 $\delta t$  – main relative permissible error limit of DE exposure time measurement that should not exceed ±5 % is defined as follows:

$$\delta t = \frac{1.1\sqrt{(\Delta t_c)^2 + (\Delta t_p)^2 + (\Delta t_o)^2}}{t}, \quad (5.7)$$

 $\Delta t_c$  - permissible error limit of the standard stop-watch;

$$\Delta t_p = 1$$
 s - error caused by a response of the user;

 $\Delta t_{o} = 1 \text{ s}$  - error caused by the source direction

towards the collimator.

5.2.7.5.9 Switch the dosimeter on and set the threshold level value of DER to 0.00  $\mu$ Sv/h, the value of DE threshold level – 0.1 mSv/h. Perform operations 5.2.7.5.6 -5.2.7.5.8 for DER  $\dot{H}_{p0}(10) = (120\pm 20) \mu$ Sv/h.

5.2.7.5.10 Switch the dosimeter on and program the threshold level value of DER to 0.00  $\mu$ Sv/h, the value of DE threshold level – 1 mSv/h. Perform operations 5.2.7.5.6 - 5.2.7.5.8 for DER  $\dot{H}_{p0}(10) = (1.2 \pm 0.2)$  mSv/h.

5.2.7.5.11 Switch the dosimeter on and set the threshold level value of DER to 0.00  $\mu$ Sv/h, the threshold level value of DE – 10 mSv/h. Then perform operations 5.2.7.5.6 - 5.2.7.5.8 for DER  $\dot{H}_{p0}(10) = (12\pm 2)$  mSv/h.

5.2.7.5.12 Switch the dosimeter on and set the threshold level value of DER to 0.00  $\mu$ Sv/h, the threshold level value of DE – 100 mSv/h. Then perform operations 5.2.7.5.6 - 5.2.7.5.8 for DER  $\dot{H}_{p0}(10) = (120 \pm 20) \text{ mSv/h}.$ 

5.2.7.5.13 Switch the dosimeter on and set the threshold level value of DER to  $0.00 \,\mu$ Sv/h, the threshold level value of DE – 8000 mSv/h.

5.2.7.5.14 Enter the "Clock" mode to adjust the current time with  $\pm 1$  min error and proceed to save the readings.

5.2.7.5.15 Enter the "DE" mode and perform operations 5.2.7.5.6 for DER  $\dot{H}_{\nu 0}(10) = (900 \pm 100) \text{ mSv/h}$ . In the period of 10 hour measurement time take DE measurement results alongside the readings in the "Clock" mode, which should differ from the stop-watch readings in not more than 5 min. Then check if the dosimeter generates audio and light signals. Perform operations 5.2.7.5.8.

5.2.7.5.16 Perform operations 5.2.7.5.6 for DER  $\dot{H}_{p0}(10) = (900 \pm 100)$  mSv/h. In the period of measurement time *t* expressed in seconds and calculated by the formula:

$$t = \left(60 + \frac{36000}{\dot{H}_{p}(10)}\right) + t_{\partial}, \qquad (5.8)$$

where  $H_p(10)$  is a numeric DER value, expressed in  $\mu$ Sv/h, take the DE measurement results. Perform operations 5.2.7.5.8. 5.2.7.5.17 Enter the "Alarm clock" mode to specify the 2 min interval with reference to real time at which the alarm clock is to be activated.

5.2.7.5.18 When the alarm clock sends audio signal, switch on the stop-watch to control the alarm clock actuation. The signal should take  $(60 \pm 2)$  s.

5.2.7.5.19 Maximum value of all received errors is the main relative permissible error limit of DE measurement. 5.2.7.5.20 The result of the dosimeter verification is considered positive if the main relative permissible error limit of DE measurement with 0.95 confidence probability does not exceed  $\pm 15$  %.

5.2.7.5.21 If the limit of main relative permissible error in DE measurement does not meet the requirements outlined in 5.2.7.5.20, the dosimeter can not be verified and should be sent for repair.

5.2.7.6 Presentation of verification results

5.2.7.6.1 Positive results of primary or periodic verification are registered as follows:

- primary verification is registered in the "CERTIFICATE OF ACCEPTANCE" section;

- periodic verification is registered in the issued Certificate of the established form.

5.2.7.6.2 If recognized inapplicable for use the dosimeter:

- should not be allowed to manufacture and use at primary verification;

- gets the certificate of inadequacy during use and after repair.

## **6 STORAGE**

6.1 The dosimeters should be stored packed, safe from mechanical damage in dry, ventilated and clean storehouses at the ambient temperature from +5 to +40 °C and relative humidity up to 80 % at +25 °C temperature. The storehouse should be free of dust, acids, gas, vapors of organic solvents, and alkali that may cause corrosion.

6.2 Maximum shelf life of the dosimeters in packing is 3 years.

## 7 SHIPPING

7.1 Packed dosimeters may be shipped at the ambient air temperature from - 30 to + 65 °C; relative humidity up to 100 % at + 25 °C temperature by railway, air, water or motor transport at any distances in conformance with the following rules:

- by railway transport in a clean box car;
- by air transport in pressurized compartments;
- by water transport in a dry hold;
- by motor transport in a closed car.

7.2 The dosimeters in shipping container should be placed and fixed in the vehicle to ensure their stable position and to avoid shocks.

7.4 Observe the inscriptions on the shipping container at loading and unloading the dosimeters.

7.5 The dosimeters must not be exposed to the influence of atmospheric precipitations during loading-unloading.

7.6 The dosimeters may be delivered in parcels in conformance with the international rules.

### 8 DISPOSAL

Disposal of the dosimeter is performed in compliance with the general rules, i.e. metals are recycled or melted, and plastic parts are dumped.

**Note.** If the dosimeter is solied by liquids or dry radionuclides, and it is impossible to completely deactivate it, it should be buried as solid radioactive waste.

## 9 WARRANTY

9.1 The manufacturer guarantees the conformity of the dosimeter to the TV V 33.2-22362867-010:2007 technical requirements provided that the customer observes the guidelines for its use, shipping and storage conditions presented in the BICT.412118.023-02 PO operating manual.

9.2 The warranty period of the dosimeter use shall terminate and be of no further effect in not less than 18 months after the date of putting it into operation, or after the end of the storage period.

9.3 The warranty period of storage of the dosimeter is 6 months after its manufacture date.

9.4 The warranty period is prolonged for the time when the dosimeter has been under warranty repair.

9.5 Warranty is invalid in case of use, shipping and storage violations, any mechanical damages, or if the warranty seals are violated. In this case the repair is performed at the user's expense.

9.6 After the warranty period terminates, the repair of the dosimeter is performed under separate contracts (as further agreed).

9.7 Warranty and post-warranty repair is done only by the manufacturer.

## **10 PACKING CERTIFICATE**

The DKG-21 M personal gamma i radiation dosimeter of BICT.412118.023-02 type with serial number is packed by the Private Enterprise "SPPE "Sparing-Vist Center" in accordance with the requirements specified in TV V 33.2-22362867-010:2007.

(position)

(signature / print full name)

(year, month, date)

#### 11 CERTIFICATE OF ACCEPTANCE

The DKG-21 M personal gamma radiation dosimeter of BICT.412118.023-02 type with \_\_\_\_\_\_serial number is manufactured to meet the technical requirements specified in TV Y 33.2-22362867-010:2007, and is accepted for use. OCD head

(signature / print full name)

Stamp here State verification officer

(signature / print full name)

Verification mark here

(year, month, date)

110

#### **12 REPAIR**

12.1 In case of failure or troubles during the warranty period of the dosimeter, the user should contact the producer enterprise by e-mail (see below) to receive the address of the nearest service center:

#### PE "SPPE "Sparing-Vist Center" Tel.: (+380 32) 242 15 15; Fax: (+380 32) 242 20 15 E-mail: sales @ ecotest.ua

## **13 ACCEPTANCE AND WARRANTY**

DKG-21 M personal gamma radiation dosimeter TY Y 33.2-22362867-010:2007

Serial number

Manufacture date


Primary calibration performed


Hereby I confirm the acceptance of the packed device applicable for use and the acceptance of the warranty terms Sales date


Salesperson signature

Stamp here

Note. If any controversy arises, the parties should act in accordance with the Art. 14 of the Law of Ukraine on the Protection of Consumer Rights

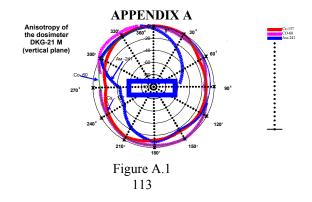

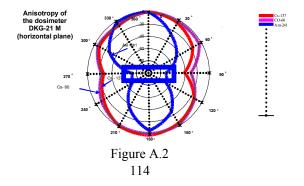

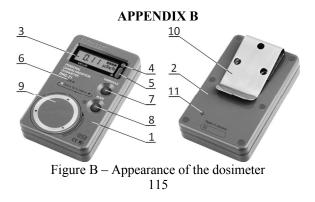

PRIMARY AND PERIODIC VERIFICATION OF KEY SPECIFICATIONS			
Verified specification		Date of measurement	
	Value according to specification	year 20	
Name		Actual value	Measured by (position, signature)
1 Main relative error limit of DER			
measurement with confidence			
probability of 0.95, % in the DER range:			
- of 1.0 μSv/h to 10 μSv/h - of 10 mSv/h to 1.0 Sv/h	±20 %		
	±15 %		

116

C-1

Date of measurement					
year 20		year 20		year 20	
Actual value	Measured by (position, signature)	Actual value	Measured by (position, signature)	Actual value	Measured by (position, signature)

PRIMARY AND PERIODIC VERIFICATION OF KEY SPECIFICATIONS			
Verified specification	Date of measurement		
	Value	у	ear 20
Name	according to specifi- cation	Actual value	Measured by (position, signature)
2 Main relative error limit of DE measurement with 0.95 confidence probability	±15 %		

C-2

Date of measurement					
year 20		year 20		year 20	
Actual value	Measured by (position, signature)	Actual value	Measured by (position, signature)	Actual value	Measured by (position, signature)